

Learn. Share. Collaborate.



ZEISS Labscope

Your microscope software to empower digital education.

zeiss.com/labscope



Seeing beyond

Your microscope software to empower digital education.

› **In Brief**

› The Advantages

› The Applications

› Technology and Details

Labscope is your easy-to-use imaging app for digital education.

Be it for the university, school or even your hobby – it's easier than ever before to snap images, record videos and measure your microscopic samples. You can easily create digital classrooms – just connect your ZEISS microscopes into a network. Explore the advantages of an interactive learning atmosphere where you can engage your students fully and enthuse them with your lessons.

No matter which device you use, you will enjoy the same look and feel: no training required. It's never been so simple and efficient until now.




▶ [Click here to view this video](#)

Simpler. More Intelligent. More Integrated.

- › In Brief
- › **The Advantages**
- › The Applications
- › Technology and Details

Enjoy Connected Microscopy

In digital education connectivity is a core feature, offering you control, remote observation and the use of mobile devices to learn, share and collaborate. With Labscope you can transform your student lab via a Wi-Fi-enabled network into a digital classroom. Labscope lets you display all the live images from all the connected microscopes at any time and from anywhere in the room. Observe how your students are doing in real-time. Project images from any microscope onto the wall or a TV screen. Discuss what students are seeing with the whole class. As you move freely around the room, 'lecture & listen' becomes 'show & see'.

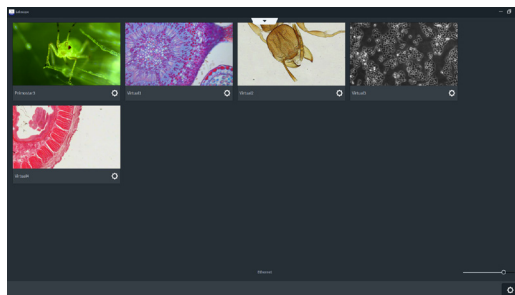

Simple. Imaging. App.

Whether you're using a traditional Windows PC with mouse and keyboard, with touchscreen, an Android device, or an iPad or iPhone we've got you covered. Labscope runs on all of these devices. Labscope leverages network technology to switch effortlessly between microscopes and get real-time thumbnails of all microscopes in the room. If you're just looking for an easy-to-use, single-microscope imaging solution, you can skip the network and connect your ZEISS microscope to any Windows computer via a USB cable.



Powerful Features with Every Touch

Get your students measuring, annotating and documenting their results by making it easy to snap images or record videos. Save your images in the ZEN-compatible .czi file format, which includes all metadata and a separate annotation layer – or select the .jpg format to save space. Use the integrated pointer to lead your students to areas of interest. Let them do their hand drawings with the drawing function. Or opt for Labscope Teacher to manage and organize your digital classroom. Labscope works easily by touch and gesture control, independently from your device.



Your Insight into the Technology Behind It


- › In Brief
- › **The Advantages**
- › The Applications
- › Technology and Details

Easy digital solutions


To connect a WiFi-enabled network that spans the whole lab or classroom, use Labscope-ready ZEISS microscopes. Such as Primostar 3 HD, Primovert HDcam, Stemi 305 cam, Axioscope 5/7/Vario and Axiolab 5 – or any other microscopes equipped with cameras like Axiocam ERc 5s, Axiocam 202 mono, Axiocam 208 color.

Document and archive your results.

And share the images in your digital network. It is full of possibilities.

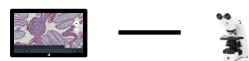


- Connect Labscope to your microscope to start with digital microscopy work.

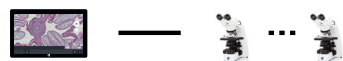


- Connect Labscope to multiple microscopes to fulfill your microscopy works on different microscopes.



- Connect your microscope to multiple tablets to do the microscopy work by different users at same time.

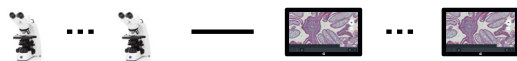


- Connect multiple microscopes to multiple devices to enable a fully connected lab or classroom.

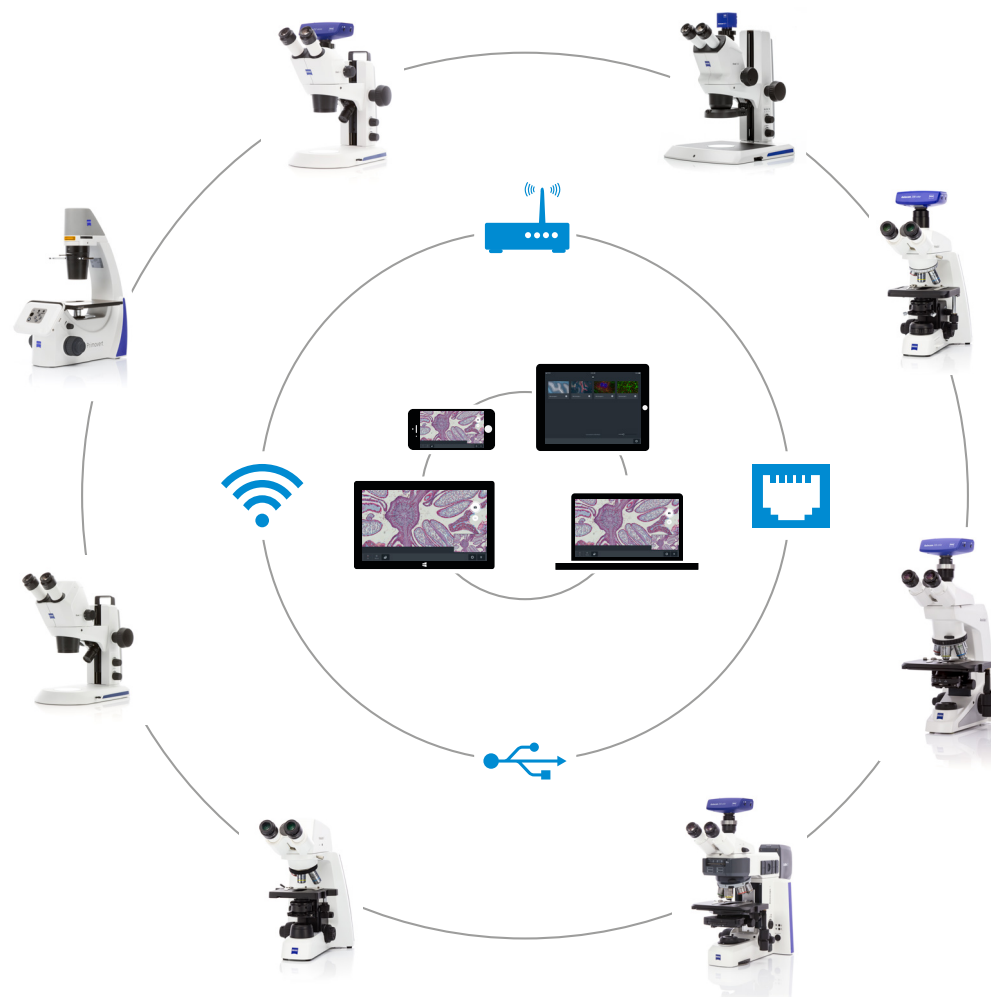

Expand Your Possibilities


- › In Brief
- › The Advantages
- › **The Applications**
- › Technology and Details

Basic functions

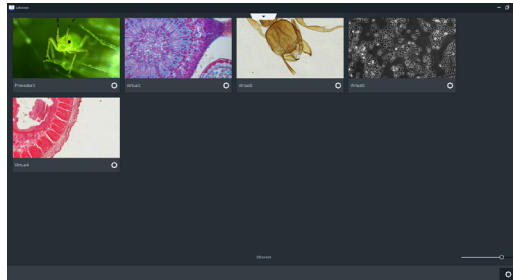


Microscope list

Get an overview of all connected microscopes. Manage and work on them easily.

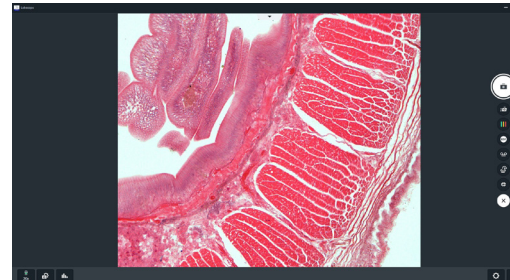


Live view and acquisition modes*

High-quality live image and multiple acquisition modes make your microscopy work more flexible.

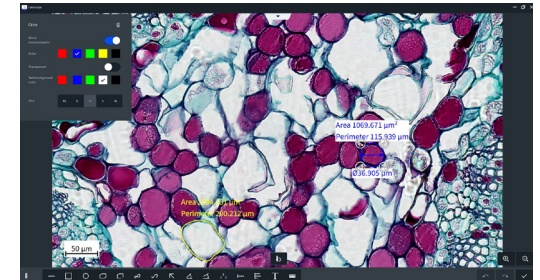


Annotations and measurements

Choose between 15 types of annotations and easy adjustments of color and size.

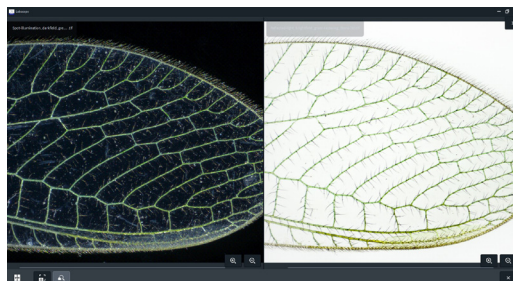


Split view

Compare your images in detail with the split view. Zoom in and zoom out, and even save the image of split view for documentation.

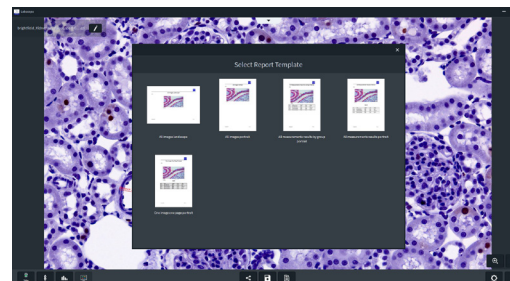

Report

Different kinds of report templates are prepared for you to quickly create reports from your microscopy work.

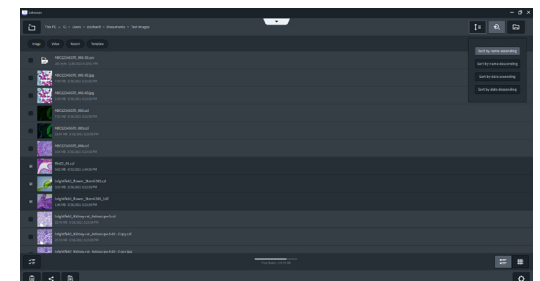

File management


Sort, search and share your images, videos and reports. Barcode scanning is also available when you want to set a file name or search a file.

* Availability of each acquisition mode is depending on cameras and microscopes.

Expand Your Possibilities


- › In Brief
- › The Advantages
- › **The Applications**
- › Technology and Details

Advanced functions

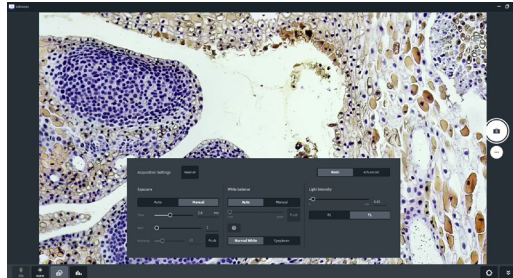

Acquisition settings

Get in-app control of your hardware to adjust exposure, white balance, light intensity and more. Focus on the specimen all the time.

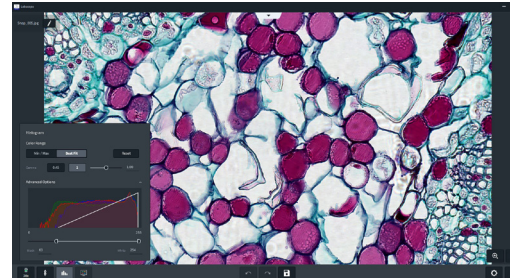


Histogram

The histogram panel is available for both live view and image view. The Min/Max and Best Fit functions optimize the display effect with just one click.




Drawing Tube

Do hand drawings of microscopy images without drawing skills. The translucent sketch solution provides an inspiring learning style.

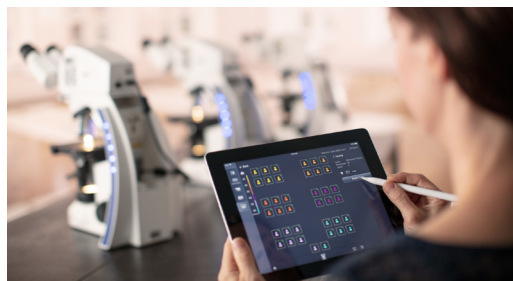

Labscope Teacher*

Organize and manage your digital classroom efficiently. The digital classroom solution from ZEISS creates an interactive learning environment where digital natives feel at home.

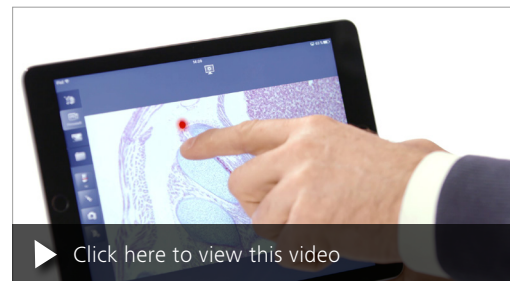

Pointer Function

Use a projector to grab student's attention and point them toward tiny structures and interesting details.

* Labscope Teacher is a licensed software, please consult with sales representative for enquiry, demo and free trial.

Your Flexible Choice of Components

- › In Brief
- › The Advantages
- › The Applications
- › **Technology and Details**

		ZEISS Labscope for iPhone	ZEISS Labscope for iPad	NEW: ZEISS Labscope for Android	ZEISS Labscope for Windows
Connections	Wireless	Wi-Fi	Wi-Fi	Wi-Fi	Wi-Fi
	Wired	–	Ethernet connection via Lightning or USB-Type-C adapters	–	USB, Ethernet
Thumbnail of microscope	Live thumbnail with adjustable size	●	●	●	●
Microscope configuration	Objective selection	●	●	●	●
	Shading correction	–	●	–	●
	Configuration lock	●	●	●	●
	Scale calibration	–	●	–	●
	QR code of configuration	●	●	●	–
Acquisition	Single snap	●	●	●	●
	Fast snap	●	●	●	●
	Video recording	●	●	●	●
	Time-lapse	–	●	–	●
	EDF	–	●	–	●
	Multi Channel*	–	○	–	○
	Fast Panorama*	–	–	–	○
Acquisition settings	Exposure, white balance, light intensity	●	●	●	●
	Adjustment of live quality, image orientation, color/grayscale switch, Gamma, denoise, sharpening, pixel correction and HDR	–	●	–	●
Live assistance	Histogram, graticule overlay, over-exposure indicator, focus indicator and split view	–	●	–	●

● functions available ○ optional – functions not available **NEW** new features in Labscope v4.01

The availability of some functions is related with types of cameras and microscopes.

* Multi Channel and Fast Panorama are licensed modules. Get more information in the module manager of Labscope and have a free trial. Please see the remarks in page 8 for supported cameras.

Your Flexible Choice of Components

- › In Brief
- › The Advantages
- › The Applications
- › **Technology and Details**

		ZEISS Labscope for iPhone	ZEISS Labscope for iPad	NEW: ZEISS Labscope for Android	ZEISS Labscope for Windows
Annotations	Line, arrow, rectangle, circle, text, count and scale bar	●	●	●	●
	Polygon, polyline, spline contour, spline, angle, disconnected angle, count, caliper and multi calipers	–	●	–	●
	Adjustment of annotations: color and size of line, text and text background	●	●	●	●
Image processing	Histogram, over-exposure indicator, split view, multi-channel comparison	–	●	–	●
	Gamma, brightness, contrast, color intensity, sharpness	●	●	–	●
Supported image types	CZI, JPEG, TIFF	●	●	●	●
File management	Different types of file views, sorting, search and sharing	●	●	●	●
	Report with multiple templates	–	●	–	●
Interfaces for external functions	Barcode scanning for configuring a file name or searching a file	–	● (iPad camera)	–	●
	Foot pedal with shortcut setting for snapping	–	● (Bluetooth foot pedal)	–	●
Theme	Dark and light theme	●	●	●	●
Global settings	General settings, default annotation settings, etc.	●	●	●	●
	Module with licensed modules	–	●	–	●
Digital classroom extensions	Pointer, drawing tube and the connection with the classroom management software Labscope Teacher*.	Pointer only	●	Pointer only	●
Extensions for iOS functions	Importing and managing Labscope files via Files App; Sharing files via AirDrop	●	●	n.a.	n.a.

● functions available ○ optional – functions not available

The availability of some functions is related with types of cameras and microscopes.

* Labscope Teacher is a licensed software. Please consult your sales representatives for an enquiry, demo and free trial.

Technical Specifications

- › In Brief
- › The Advantages
- › The Applications
- › **Technology and Details**

Application Information:	
Seller	Carl Zeiss AG
Category	Education, Documentation, Microphotography, Laboratory, Research
Compatible ZEISS microscopes	Primostar 3 HD, Primotech (iOS and Windows only), Primovert HDcam, Stemi 305 cam, Axioscope 5/7/Vario and Axiolab 5
Compatible ZEISS cameras	Axiocam ERc 5s, Axiocam 202 mono, Axiocam 208 color, Axiocam 105 color (Windows only) and Axiocam 305 mono/color (Windows only)
Languages	English, Czech, French, German, Italian, Japanese, Korean, Polish, Portuguese, Russian, Simplified Chinese, Spanish




System requirements	ZEISS Labscope for iPhone	ZEISS Labscope for iPad	NEW: ZEISS Labscope for Android	ZEISS Labscope for Windows
Operating system	iOS 13 or later	iOS 13 or later	Android 9.0 or later	Windows 10 (64-bit)
Minimum hardware	iPhone 6s/plus iPhone SE (2016)	iPad Air 2 iPad 5 iPad Mini 4	RAM: 4 GB	CPU: i3 (6 th generation) dual-core @ 2.5 GHz RAM: 4 GB
Recommended hardware	iPhone 7 or later	iPad Air 3 or later iPad 6 or later iPad Mini 5 or later iPad Pro 1 or later	RAM: 6 GB or above	CPU: i5 (7 th generation) quad-core @ 3.0 GHz or above RAM: 8 GB or above

System requirements	ZEISS Labscope Module Fast Panorama
Operating system	Windows 10 (64-bit)
Minimum hardware	CPU: i5 (8 th generation) quad-core @ 3.0 GHz RAM: 8 GB or above
Recommended hardware	CPU: i7 (9 th generation) hexa-core @ 4.0 GHz or above RAM: 16 GB or above

Supported camera types for Multi Channel module: Axiocam 105 color, Axiocam 202 mono, Axiocam 208 color and Axiocam 305 mono/color.

Supported camera types for Fast Panorama module: Axiocam 305 mono/color.

Not all performance specifications of the used cameras can be supplied in the Labscope environment due to application specific restrictions.

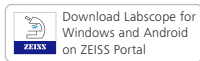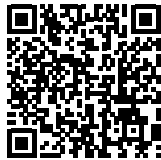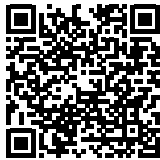

*Download Labscope for
Windows and Android*

*Download Labscope
for Android*

*Download Labscope
for iOS*

Note:

For service & support visit: [ZEISS Labscope Forum](#).

For ideas of product development: [ZEISS Portal My Voice](#).



Carl Zeiss Microscopy GmbH
07745 Jena, Germany
microscopy@zeiss.com
www.zeiss.com/labscope